

Insta-TEST[®]

25 Water Test Strips
Low Range Peroxide

Just DIP & READ!



LaMotte
 Chestertown, MD 21620 USA
 www.lamotte.com



Immerse test strip. Remove with pad face up.
DO NOT SHAKE OFF EXCESS WATER.
 Immediately compare to color chart.

Peroxide ppm	Code 2984LR
0	
1	
3	
10	
30	
50	

- Keep wet fingers out of the vial.
- For best results, immerse strip to a depth of 12" - 18".
- Close vial tightly after removing strip.
- Store in a cool, dry place.

