




Peracetic Acid
ppm

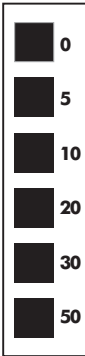

TO USE:

Using a cup-size sample, immerse test strip and remove with **pads face up. DO NOT SHAKE OFF EXCESS WATER.** Wait 10 seconds and compare to color chart immediately. Discard strip after 30 seconds.



TIPS

- Keep wet fingers out of the vial.
- Close vial tightly after removing strip.
- Store in a cool, dry place.

LaMotte

Chestertown, MD 21620
<http://www.lamotte.com/insta>

made in
USA★

Code 3000-LR