## Rapid Test for determination of protein residues

### Contents:

- 1 Aluminium container with 60 test strips
- 1 Bottle INDIPRO-1
- 1 Bottle INDIPRO-2

INDIPRO is a rapid test for determination of protein residues on working surfaces and tools having contact with foodstuffs.

### Safety precautions:

Reagent INDIPRO-1 contains ethanol 35–55%, Reagent INDIPRO-2 contains ethanol 20–35% CAS 64-17-5. **WARNING.** Flammable liquid and vapor. Keep away from heat/sparks/open flames/hot surfaces. – No smoking. Keep container tightly closed. Store in a well-ventilated place. Keep cool. With reduced labelling acc. CLP Annex I chapter 1.5.2.1. For further information, please ask for safety data sheet (*see [www.mn-net.com/MSDS](http://www.mn-net.com/MSDS)*).

### General indications:

Remove only as many test strips as are required. Close the container immediately after removing a strip. Do not touch the test field.

### Instructions for use:

1. Add 1 drop of reagent INDIPRO-1 on the test field of the test strip.
2. Draw test strip with the moistened test field approximately 10 cm under low pressure over the surface to be tested.
3. Add 1 drop of reagent INDIPRO-2 on the test field. A colour change from yellow to green indicates the presence of protein residues. The surface has to be cleaned.

### Limit of sensitivity:

50 µg BSA (bovine serum albumin) gives a clear green colouration

### Interference:

Residues of desinfectants can cause also a green colouration of the test field.

### Disposal:

Dispose of used strips at an authorised hazardous waste collection point.

### Storage:

Avoid exposing the strips to sunlight and moisture. Keep container cool and dry (storage temperature not above +30 °C).

If correctly stored, the test strips may be used until the use-by-date printed on the packaging.

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